

---

Poster

## [P25-9] P25-9: Oncologic drugs (1)

Chair: Ryuji Ikeda, Japan

Mon. Sep 25, 2017 12:30 PM - 1:30 PM Annex Hall (1F)

---

(Mon. Sep 25, 2017 12:30 PM - 1:30 PM Annex Hall )

## [P25-9-9] UPLC-MS/MS analysis of plasma busulfan for clinical research

Gareth Hammond<sup>1</sup>, Stephen Balloch<sup>2</sup> (1.Waters Corporation, 2.Waters Corporation)

Keywords: Busulfan, Plasma, Mass Spectrometry, LC-MS/MS

### Background

Busulfan is a bifunctional alkylating agent whose bioavailability varies greatly between individuals due to factors such as age, underlying diseases and drug-drug interactions. An accurate, sensitive and specific analytical method may play a role in assessing the pharmacokinetic and pharmacodynamic effects of busulfan administration in clinical research.

### Methods

Samples (50L) were deproteinised with busulfan-<sup>2</sup>H<sub>8</sub> internal standard in methanol. Elution was achieved within 3 minutes using a HSS-T3 C18 UPLC column (2.1x50mm, 1.8m) on the Waters ACQUITY UPLC<sup>®</sup> I-Class with a water/methanol/ammonium acetate/formic acid gradient. Ammonium adducts of busulfan were analyzed using electrospray ionization in positive mode with multiple reaction monitoring using the Waters XEVO<sup>®</sup> TQD mass spectrometer.

Matrix-matched calibrators (0.025-5 g/mL) and quality control samples (0.05, 0.75, 1.5 and 3.5 g/mL) were prepared. Due to known instability of busulfan in plasma, aliquots of in-house calibrators and quality control samples were stored frozen and thawed prior to use each time the method was used.

### Results

Analytical sensitivity was calculated to be 0.020 g/mL (n=10 extractions, five occasions, 16.0% CV). Linearity was demonstrated over the concentration range 0.0175-6.51 g/mL and system carryover was negligible in samples 10 g/mL.

Precision studies (n=5, five occasions) demonstrated repeatability and total precision 7.3%. A comparison was made by analysing anonymized plasma samples (n=40, range 0.20-2.28 g/mL) against an independent UPLC-MS/MS method. An ordinary linear fit comparison of r=0.998, a mean bias of 5.3% from Altman-Bland analysis and a Deming equation of  $y = 1.01x + 0.04$  were obtained.

The mean recovery for busulfan pooled plasma samples (n=3, 0.05 and 3.5 g/mL) was between 85.1-106.1% in the presence of high concentrations of endogenous compounds albumin, bilirubin, cholesterol, triglycerides and uric acid and exogenous Intralipid<sup>®</sup>. The mean recovery was between 93.2-103.0% in the presence of acetaminophen, fluconazole, ketoconazole, itraconazole, phenytoin, posaconazole and voriconazole.

Negligible matrix effects were observed at low (0.05 g/mL) and high (3.5 g/mL) concentrations as indicated by respective mean internal standard adjusted matrix factors of 0.99 (range 0.97-1.02) and 0.98 (range 0.95-1.00).

### Conclusions

The clinical research method developed for the quantification of plasma busulfan demonstrated good linearity, analytical sensitivity and precision with negligible matrix effects.

For Research Use Only, Not for use in diagnostic procedures.